

TransScript® IV Reverse Transcriptase [M-MLV, RNase H-] (Anti Inhibition High Temperature RT)

Please read the data sheet carefully prior to use.

Cat. No. AW101

Version No. Version 1.0

Storage: at -18°C or below for two years

Concentration: 200 units/μl

Description

TransScript® IV Reverse Transcriptase is an M-MLV reverse transcriptase with high thermostability expressed and purified from *E. coli* after genetic modification. It enables a board range reaction temperature (42°C-65°C) and good resistance to inhibitors, with ultra-high thermal stability, and deficient RNase H activity. Its high reaction temperature is beneficial to open RNA secondary structure when synthesizing the first strand cDNA, with the optimal reaction temperature of 50°C. It also has the characteristics of **good resistance to inhibitors**, **fast reaction speed**, high sensitivity, high specificity, high yield (more full-length cDNA), high thermostability, and long half-life.

Features

- Wide range of applications: severely degraded RNA can be used as a template for reverse transcription.
- Agnosticto inhibitors remaining in RNA.
- Fast reaction speed: only **10** minutes for reverse transcription.
- Strong synthesis ability: synthesis length up to 20 kb.
- High thermal stability: reaction temperature 42°C-65°C.
- High sensitivity: high detection rate for trace RNA.

Applications

- High-copy number and low-copy number gene detection.
- High GC or complex secondary structure RNA template.
- For difficult-to-handle RNA sample: degraded RNA template and RNA template containing RT enzyme inhibitor.
- cDNA library construction, primer extension, 3' and 5' RACE.

Kit Contents

Component	AW101-02
TransScript® IV Reverse Transcriptase	10000 units
10×TS IV RT Buffer	100 μl
Anchored Oligo (dT) ₂₀ Primer (0.5 μg/μl)	50 μl

Prior to use, please centrifuge each component briefly.

First-strand cDNA synthesis

1. Reaction components

Component	Volume
Total RNA/mRNA	0.1 ng-5 μg/ 10 pg-500 ng
Anchored Oligo(dT) ₂₀ Primer (0.5 μg/μl)	1 μl
or Random Primer (0.1 μg/μl)	1 μl
or GSP	2 pmol
10 mM dNTPs	1 μl
10×TS IV RT Buffer	2 μl
Ribonuclease Inhibitor	0.5 μl
TransScript® IV Reverse Transcriptase	1 μl
RNase-free Water	Variable
Total volume	20 μl



2. Mix well gently

- If use Anchored Oligo (dT)₂₀ or GSP, incubate at 50°C for 10 minutes.
- If use Random Primer (N9), incubate at 25°C for 10 minutes, and then incubate at 50°C for 10 minutes.
- For GC-rich or complex secondary structure RNA template, it is suggested to increase the reaction temperature to some extent (≤65°C).

3. Incubate at 85°C for 5 seconds to inactivate *TransScript*[®] IV Reverse Transcriptase.

Recommended qPCR system and conditions (taking 20 µl reaction system as an example)

Component	Volume	Final Concentration
Template	Variable	As required
Forward Primer (10 µM)	0.4 µl	0.2 µM
Reverse Primer (10 µM)	0.4 µl	0.2 µM
2× <i>PerfectStart</i> [®] Green PCR SuperMix	10 µl	1×
Passive Reference Dye (50×) (optional)	0.4 µl	1×
Nuclease-free Water	Variable	-
Total volume	20 µl	-

qPCR (Three-step)

94°C 30 sec
 94°C 5 sec
 50-60°C 15 sec ★
 72°C 10 sec ★

40-45 cycles

Dissociation Stage

qPCR (Two-step)

94°C 30 sec
 94°C 5 sec
 60°C 30 sec ★

40-45 cycles

Dissociation Stage

For the ABI qPCR instrument, we suggest using the following read time (Fluorescent signals can be collected during the annealing or extension stage for three-step qPCR):

- ★ For ABI Prism7700/7900, set the read time to 30 seconds.
- ★ For ABI Prism7000/7300, set the read time to 31 seconds.
- ★ For ABI Prism7500, set the read time to 34 seconds.
- ★ For ABI ViiA7, set the read time is at least 19 seconds.

Three-step qPCR is more suitable for higher amplification efficiency assay.

Two-step qPCR is more suitable for higher specificity assay.

Recommended PCR system and conditions (taking 50 µl reaction system as an example)

Component	Volume	Final Concentration
Template	Variable	as required
Forward Primer (10 µM)	1 µl	0.2 µM
Reverse Primer (10 µM)	1 µl	0.2 µM
2× <i>TransTaq</i> [®] HiFi PCR SuperMix II	25 µl	1×
Nuclease-free Water	Variable	-
Total volume	50 µl	-

PCR

94°C 2-5 min
 94°C 30 sec
 50-60°C 30 sec
 72°C 1-2 kb/min
 72°C 5-10 min

35-40 cycles



Notes

- Avoid RNase contamination.
- For complex RNA templates, or to obtain higher synthesis efficiency, it is recommended to mix the RNA template and primers with RNase-free Water well, incubate at 65°C for 5 minutes, and put on ice for 2 minutes before adding other reaction components.
- Mixing all the reaction components in one step can complete most reverse transcription reactions. For complex RNA templates, or to obtain higher synthesis efficiency, it is recommended to add thermal incubation steps for the template and primers according to the manual.

For research use only, not for clinical diagnosis.

Version number: V1-202008

Service telephone +86-10-57815020

Service email custserv@transgenbiotech.com

