

Low Speed Centrifuge

Packing list

Main unit ×1, plastic rotor and supplied adaptor × 2, Simple user manual

A8-2P

A4-PCR8P



2

SA05P2


Installation Instructions

1. Unpack and place the centrifuge on a solid and flat surface (Laboratory bench space).
2. Ensure good working environment and ambient temperature of 15°C-25°C.
3. Protect the centrifuge from direct sunlight or temperature of higher than 30°C.
4. Place the unit at >10cm distance from other objects to allow good ventilation& performance.
5. Please keep the centrifuge away from heat or water (liquid) and chemicals.

Rotor installation

1. Ensure the centrifuge chamber is clean and dry.
2. Ensure there is no sign of corrosion or crack or damage on the rotor.
3. Place the rotor in position on the rotor base or motor shaft.
4. Press the rotor inward on the motor shaft to secure tightly.
5. Please use your both hands to load the rotor.



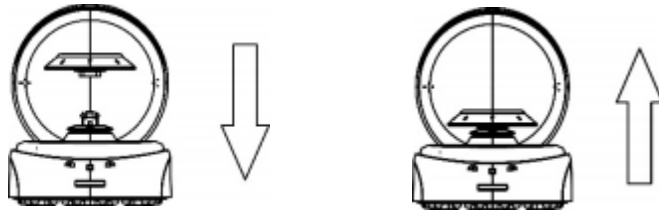

Rotor installation

1. Switch power ON located at the back of the unit.
2. Open the lid and load the centrifuge tubes symmetrically on to the rotor.
3. Close the centrifuge lid to start centrifugation process.
4. Switch power 'OFF' or open the lid and the rotor will stop gradually.
5. Open the lid and collect the samples.

202502

