

Trans5 α Chemically Competent Cell

Please read the manual carefully before use

Cat. No. CD201

Version No. Version 1.1

Storage: at -70°C or below for one year. Do not store in liquid nitrogen.

Description

Trans5 α Chemically Competent Cell is specifically designed for chemical transformation of DNA. It permits a transformation efficiency of over 10⁸ cfu/ μ g DNA (tested by pUC19 plasmid DNA).

Genotype

F ϕ 80d *lacZ* Δ M15 Δ (*lacZYA-argF*) U169 *end* A1 *recA1* *hsdR17* (*r_k⁻*; *m_k⁺*) *supE44* λ -*thi-1* *gyrA96* *relA1* *phoA*

Features

- High transformation efficiency: >10⁸ cfu/ μ g (pUC19 DNA).
- Reduced recombination of cloned DNA.
- Blue/white selection.

Procedures

- Equilibrate a water bath to 42°C.
- Warm a vial of SOC medium or LB medium to room temperature. Warm selective plates at 37°C for 30 minutes.
- Thaw a vial of 100 μ l of Trans5 α Chemically Competent Cell on ice, aliquot 50 μ l of the cells into a prechilled 1.5 ml tube, add target DNA (1 to 5 μ l) into the tube. Do not mix by pipetting up and down. Incubate the cells on ice for 30 minutes.
- Heat-shock the cells for 45 seconds at 42°C without shaking. Immediately transfer the tube to ice. Incubate on ice for 2 minutes without shaking.
- Add 500 μ l of prewarmed SOC medium or LB medium (without antibiotic) into the tube, mix well and shake at 37°C for 1 hour at 200 rpm for cell recovery and for the expression of antibiotic resistance.
- Spread 20 to 200 μ l from each transformation vial on a prewarmed selective plate. The remaining can be stored at 2-8°C and plated the next day if needed.
- Invert the plate and incubate at 37°C overnight.
- Select colonies and analyze by restriction enzyme digestion, PCR, or sequencing.

Notes

- Higher efficiency transformation can be achieved by transforming cells immediately following thawing.
- Avoid repeated thawing.
- Gentle handling is required for the entire procedure.

For research use only, not for clinical diagnosis.

Version number: V1.1-202512

Service telephone +86-10-57815020

Service email custserv@transgenbiotech.com

